



BELCO GLASS, INC. USA
340 Edrudo Road
Vineland, NJ 08360

BELCO ROLL-IN INCUBATORS

:

MAINTENANCE:

NOTE: Disconnect the power cord from the power source before performing any service or maintenance on this unit.

- 1) **Cleaning** Clean the incubator interior & remove & clean shelves on a regular basis. Use a disinfectant that is suitable for your application. A thorough periodic cleaning is strongly recommended. Use care when cleaning the door gasket to prevent damage which could impair the positive door seal.

Bellco recommends the following cleaning protocol:

- 2) Wash the cabinet exterior with a mild soap and water solution. Rinse with clear water, then wipe dry. Sanitizing solution such as Roccal (National Laboratories) can be used to disinfect the interior.

Caution: Never clean the unit with alcohol or flammable cleaners with the unit connected to the electrical supply. Always disconnect the unit from the electrical service when cleaning & assure all volatile or flammable cleaners are evaporated & dry before attaching the unit to the power supply.

Storage:

To prepare the unit for storage, remove all shelves & shelf clips & disconnect the power supply. Be certain that the chamber is clean & dry & the door is in the close position. For lifting / handling & proper transport procedures – see below

Location:

In locating the incubator, consider all possible conditions that might affect its performance. Do not place near extreme heat (steam radiators, stoves, ovens, etc.). Avoid direct sun rays and rapidly moving air currents.

Moving:

This unit should be moved upright on its original shipping pallet using a forklift or pallet jack. Care should be taken not to tilt the unit while on pallet. Slide unit into place.

WARNING: Do not stand behind unit when pushing from pallet! Do not tip unit in excess of 20 degrees! Push only on edges of door! Do not push on glass!

No maintenance is required on the electrical components. If electrical or temperature problems occur, please see the trouble shooting guide in the manual or contact Bellco Service Dept @ 1-800-257-7043

Helpful tips:

In the event the incubator does not operate properly, check the following before calling for Bellco's service department:

- 1. Is electrical power reaching the unit? (Check for blown fuse.)**
- 2. Is service cord connected to outlet?**
- 3. Is door closed tightly?**