



**BELCO GLASS, INC. USA**  
340 Edrudo Road  
Vineland, NJ 08360

## **Cleaning Silicone & Rubber Components:**

Silicone & rubber components may absorb cleaners & disinfectant solutions which may leak out during use. It is best not to use these cleaners or solutions to clean these products. After use, place silicone & rubber components in distilled water. Let soak. Rinse exhaustively with tap water. Be sure tubing is completely flushed. Place items in a container of distilled water & put through a wet autoclave cycle or boil. Rinse items in Distilled or deionized water. Inspect. Stained or soiled items should be discarded.